sartorius

**Installation Instructions | Installationsanleitung
Notice d'installation | Manuale d'installazione
Instrucciones de instalación**

**Sartorius
YCC01-USBM2,
YC012 (USB-SBI), YC013 (USB-BPI),
YCRI-CC13**

RS-232/USB Interface Cable
Verbindungskabel RS232/USB
Câble de connexion RS232/USB
Cavo di collegamento RS232/USB
Cable de conexión RS232/USB

98647-003-67

English page 3

In cases involving questions of interpretation, the German-language version shall prevail.

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Contents

Contents	3
Intended Use	3
Compatibility	4
Warnings and Safety Precautions	4
Installation	5
Initial Connection	6
Notes on Installation	7
Declaration of Conformity	33

Intended Use

This USB interface device cable is a solution that enables connection of a balance or scale to a computer via the USB (universal serial bus) port. The USB interface is implemented as a virtual serial interface (virtual COM port). For the application program running on the computer, there is no difference between a serial interface physically installed in the PC and one enabled in the form of a virtual interface. The interface offers the same capabilities in either case.

Compatibility

- YCC01-USBM2: Connect a Sartorius weighing instrument with 25-pin data interface
- YCO12 (USB-SBI), YCO13 (USB-BPI): Connect a Sartorius PMA series scale
- YCRI-CC13: Connect a Sartorius laboratory refractometer

System Requirements

- Computer (PC) running Windows® 98SE, Windows® ME, Windows® 2000, Windows® XP, Windows® Vista, Windows® Server 2008, Windows® 7 or Windows® 8/8.1 operating systems
- Available USB port

Warnings and Safety Precautions

Installation

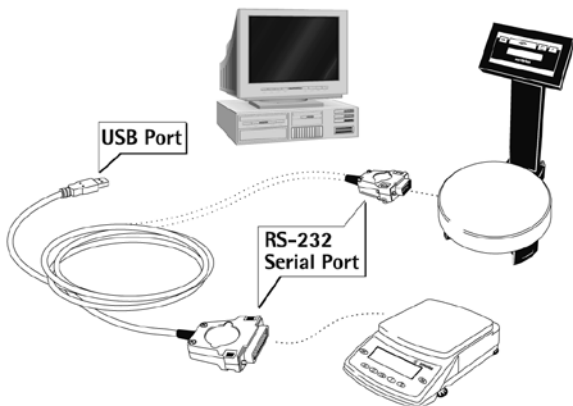
- ⚠ Be sure to turn off the power to the equipment on either side of the cable before connecting or disconnecting this cable.
- Securely fasten the D-SUB male connector to the balance or scale.

Safety

- ⚠ Use the cable only indoors and only in such a way that penetration by foreign objects or liquids is prevented.
- ⚠ Use the USB interface connector only outside hazardous areas.

Installation

- Disconnect the balance scale power supply:
Unplug the AC adapter from the wall socket (mains).

- Connect the cable to the balance scale and to the USB port on the computer.
 - Plug in the balance scale and switch it on.
- > Windows® detects the cable connected to the USB port. The first time the cable is detected, Windows® automatically runs the Wizard for adding/removing hardware.

Initial Connection

Loading the Software Drivers

Download the VCP driver online: <http://www.ftdichip.com/FTDrivers.htm>

Once you have downloaded the software, set up the virtual port on the computer.

Installation Guides

The installation guides for the drivers can be found here: <http://www.ftdichip.com/Documents/InstallGuides.htm>
Installing the Drivers for the USB Interface

Installing the Software Drivers for the USB Interface

1. Start the installation process for the downloaded program.
2. Each version of Windows® may have slightly different procedures for loading the drivers; however all versions have a wizard.

Windows® usually adds the virtual interface in the position following your highest-numbered COM port. Insert paragraph spacing as per source. For example, on a PC with 4 COM ports, the virtual interface is added as COM5 (shown in the Device Manager).

Notes on Installation

Changing the Port Number

If you use the USB interface with a program that limits the number of COM port designations (e.g., only COM1 through COM4), you may have to assign one of these port numbers to the new virtual interface.

- Set up is as follows in Windows:
Select “Start” → “Settings” → “Control Panel”
- Select “System”
- Select “Hardware” Insert arrow “Device Manager”
- Expand “Ports
- Double-click on “USB Serial Port
- Select “Port Settings” Insert arrow “Advanced”

Changing the Latency Timer

In the dialog opened above for changing the port number, you can set the latency timer to 1 msec for faster data communication.

Plug & Play Mode with Autoprint (SBI)

To use the autoprint function, you need to deactivate the plug & play mode. This setting is configured in the same dialog as that described above for changing the port number.

Using Different USB Ports on a Single PC

Driver installation must be performed each time the cable is connected to a different USB port. It is a good idea to use the same port every time, if possible.

Deinstalling the Driver

- Select "Start" → "Settings" → "Control Panel"
- Select "Software"
- Select "FTD1USB Serial Converter Driver"
- Click on "Add/Remove"
- > The software is deinstalled.

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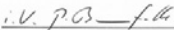
to which this certification relates is in conformity with the following standard(s) or other normative document(s) (see page 2) pursuant to the provisions of the "Directive 94/9/EC of the European Parliament and the Council of 23 March 1994 on the approximation of the laws of the Member States concerning equipment and protective systems intended for use in potentially explosive atmospheres". This product is labelled as follows:

II 3G Ex nA II T4 X

Sartorius Weighing Technology GmbH
Goettingen, 2011-10-13

Dr. Reinhard Baumfalk
Vice President R&D

Dr. Dieter Klausgrete
Leitung International Certification Management
Head of International Certification Management

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This declaration certifies conformity with the above mentioned EC Directives, but does not guarantee product attributes. Unauthorised product modifications make this declaration invalid. The safety information in the associated product documentation must be observed.

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Compliance with the Essential Health and Safety Requirements has been assured by compliance with:

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Explosive atmospheres - Part 0: Equipment - General requirements

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Explosive atmospheres - Part 15: Equipment protection by type of protection „n“

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Safety instructions:

1. *Connect or disconnect the cable only when the equipment is not energized on either side of this cable.*
2. *Securely fasten the D-SUB male connector to the PMA750...Y... by tightening the screws.*
3. *Install the cable only below buildings so that penetration of liquids or ingress of foreign bodies (such as dust or particles) is prevented.*
4. *Use the USB connector only outside the hazardous location/area.*

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This Declaration of Conformity has been drawn up on the basis of the in-house test report SAG.05.ATEX.003 written by Sartorius AG of Goettingen, Germany.

* * * * *

Sartorius Lab Instruments GmbH & Co. KG
Weender Landstrasse 94–108
37075 Goettingen, Germany

Phone +49.551.308.0
Fax +49.551.308.3289
www.sartorius.com

Copyright by Sartorius, Goettingen, Germany.
All rights reserved. No part of this publication may be
reprinted or translated in any form or by any means without
the prior written permission of Sartorius.

The status of the information, specifications and
illustrations in this manual is indicated by the date given
below. Sartorius reserves the right to make changes to
the technology, features, specifications and design of the
equipment without notice.

Status:

December 2013, Sartorius Lab Instruments GmbH & Co. KG,
Goettingen, Germany

Printed in the EU on paper bleached
without chlorine. | KT
W_YCC01-USBM2_YC012
Publication No.: WYC6029-p131206

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Diese Erklärung wird erstellt auf Basis des internen Prüfberichts SAG.05.ATEX.003 der Sartorius AG Göttingen.